


**Instruction Sheet
for the PASCO
Model EM-8648**

MAGNET SET OF TWO

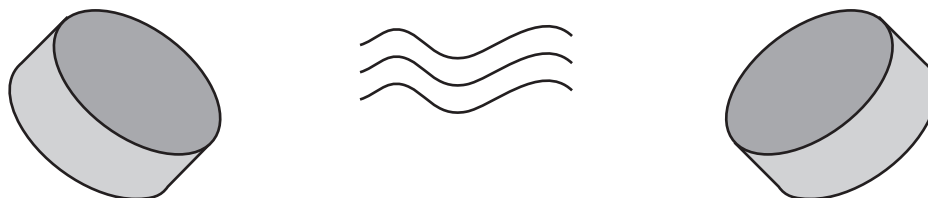

CAUTION

The PASCO Magnet Set of Two are extremely strong. When opposite poles are brought close to each other they will accelerate rapidly and easily pinch fingers. They can also erase computer discs and distort computer monitors and TV sets. We suggest you keep these magnets away from computers and your data storage.



Storing magnets pole piece to pole piece makes them difficult to separate.

Figure 2

Introduction

The PASCO Magnet Set of Two consists of two neodymium magnets suitable for a variety of experiments involving magnetism.

Storing Magnets

When storing magnets place them in a side by side orientation (see Figure 1).

It does not damage the magnets to store them pole piece to pole piece, but it makes them quite difficult to separate (see Figure 2).



Preferable to store magnets side to side.

Figure 1

Limited Warranty

PASCO scientific warrants this product to be free from defects in materials and workmanship for a period of one year from the date of shipment to the customer. PASCO will repair or replace, at its option, any part of the product which is deemed to be defective in material or workmanship. This warranty does not cover damage to the product caused by abuse or improper use. Determination of whether a product failure is the result of a manufacturing defect or improper use by the customer shall be made solely by PASCO scientific. Responsibility for the return of equipment for warranty repair belongs to the customer. Equipment must be properly packed to prevent damage and shipped postage or freight prepaid. (Damage caused by improper packing of the equipment for return shipment will not be covered by the warranty.) Shipping costs for returning the equipment, after repair, will be paid by PASCO scientific.



PASCO scientific

10101 Foothills Blvd.
Phone (916) 786-3800

P. O. Box 619011
FAX (916) 786-8905

Roseville, CA 95661-9011 USA
TWX 910-383-2040